



800-293-3003

www.emssales.net

CLEARING SOLUTION FORMULA

****CAUTION** HAZARDOUS MATERIALS**

(SPECIAL CARE SHOULD BE EXERCISED WHEN USING THESE CHEMICALS)

PLEASE REFER TO MSDS ON CHEMICALS MENTIONED BELOW FOR SAFE HANDLING

1. Add 5 milliliters of Dimethyl Phthalate
2. Add 5 milliliters of Diethyl Oxalate
3. Add one sheet (+/-0.5 gram) Filter Media 0.8 μ MCE
4. Stir mixture until Filter Media has completely dissolved
5. Before each use, check solution on a prepared blank slide to ensure that contamination or crystal formation has not occurred

****NOTE** Refrigeration of the combined chemicals greatly extends the expected shelf life. Each sheet has been cut to weigh approximately 0.5 grams (0.45-0.55)**